Pittcon 2008
59th Pittsburgh Conference on Analytical Chemistry and Applied Spectroscopy

March 1st - March 7th, 2008 Ernest N. Morial Convention Center New Orleans, Louisiana, USA

Solid Phase Extraction using Molecular Recognition Technology for Highly Selective Platinum Group Metals Separations

Session 1290 - Sample Preparation: General Day and Time: Tuesday, March 04, 2008, Morning

Molecular Recognition Technology – Solid Phase Extraction (MRT-SPE)

E-mail to: furusho@gls.co.jp

Abstract

Platinum Group Metals (PGMs) have scarcity value, and are broadly used as purification catalysts in fuel cells and diesel engine oils in recent years. The demand of PGMs for these and other applications is increasing.

However, there is concern that, in the future, the production of PGMs may not keep up with this increasing demand. As a result, recycling of PGMs from spent fuel catalysts or from base metal by-product refining is becoming important.

For the separation of PGMs for analysis, traditional methods use sedimentation, liquid-liquid-extraction, or ion-exchange extraction. Such techniques require many extraction steps and are labor-intensive.

In this study, an efficient way to separate PGMs was demonstrated using solid phase extraction (SPE) with molecular recognition technology (MRT). For the separation resin, AnaLig® PM-01, 05, 07 and 08 were used, and the adsorption behavior of PGMs to AnaLig® and the elution performance after rinse steps were examined.

Non-metal needle modified ASPEC XLi, which is an automatic solid phase extraction pretreatment method, was used for the metal analysis. After SPE extraction, the metal extracts were injected into inductivity coupled plasma optical emission spectroscopy (ICP-OES), after which the recovery rate of PGM from the AnaLig® was calculated. Finally, the selectivity of this method versus interfering metals was examined. AnaLig® PM-08 was found to be highly selective for PGMs in acid solution.

Background and Objective










Purification of Platinum Group Metals (PGMs) by SPE

Re-Cycling demand increase about PGM

Current ICP-OES technique is poor for this topics

Develop a useful method for PGM analysis using MRT-SPE-ICP-OES

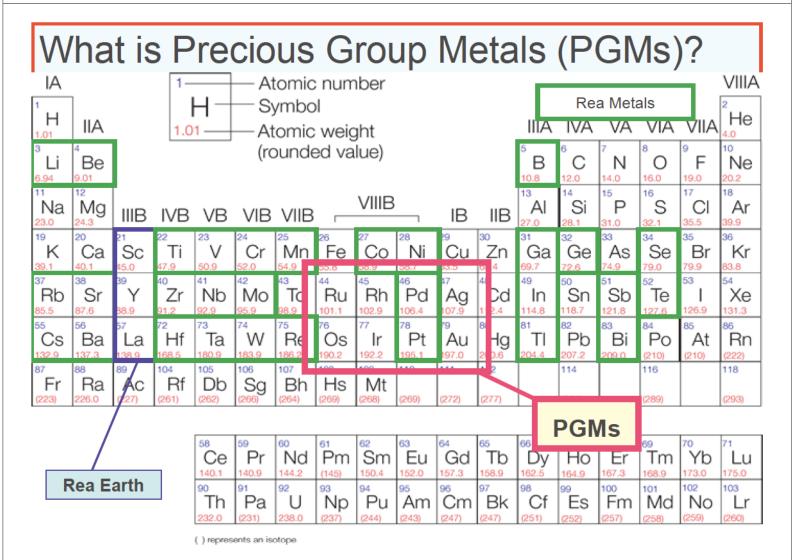

MRT-SPE Method for PGMs from High Matrix

Typical current Technique

Open column with ion Ex resin for separation of PGMs

Anion Ex resin is not good for high selective separation from matrix

Not enough clean up for determination of PMGs by ICP-OES

Spectral Interference Physical Interference Chemical Interference Present method


Separate PGMs by MRT-SPE

SPE cartridge is much easier handling than open column


MRT-SPE can separate target PGM element from main metal matrix


Reduce each interference

MRT-SPE can separate PGMs

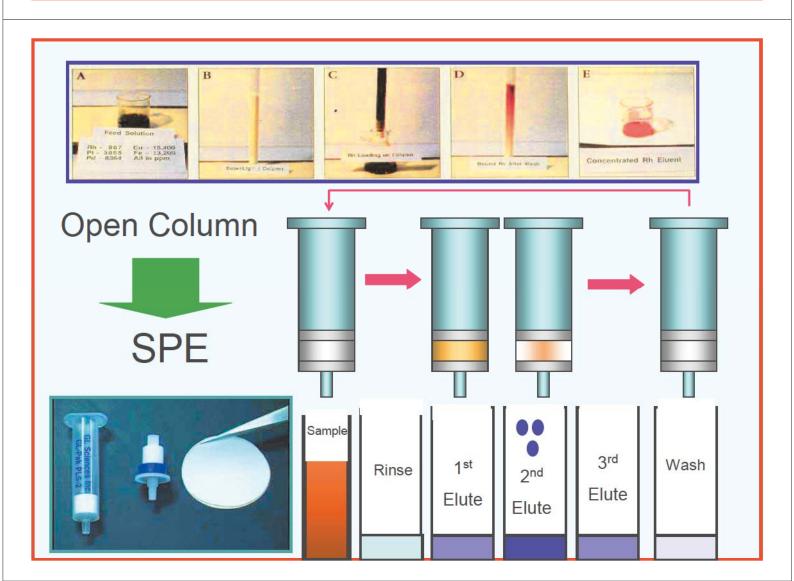


SPE Manifold System for Metal Separation

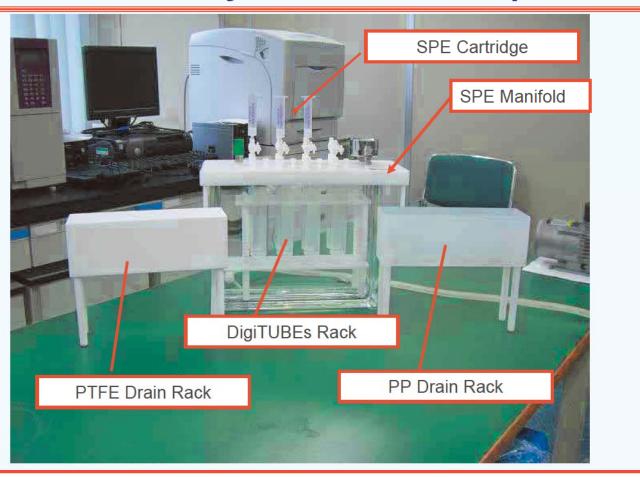

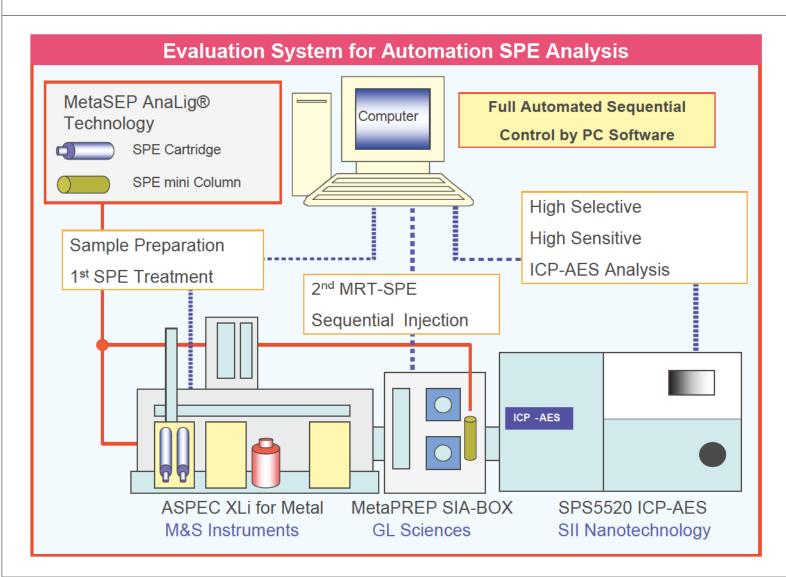



Method Development using Automated SPE system

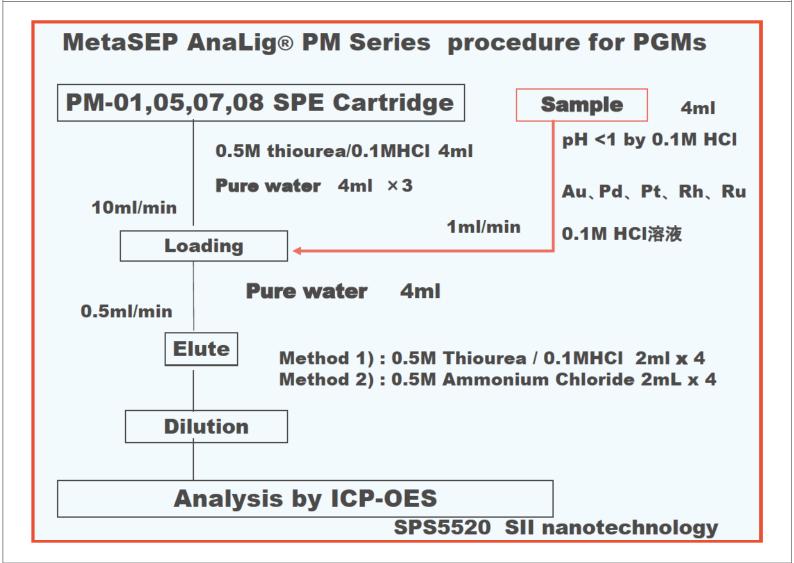

Recovery Test of MetaSEP AnaLig PM Series for PGMs

Cartridge size: 500mg/3mL

	Au		Pd		Pt		Rh		Ru		lr	
	BT	EL	ВТ	EL	ВТ	EL	ВТ	EL	ВТ	EL	ВТ	EL
PM-01	0.0%	99.5%	0.4%	88.0%	1.4%	76.9%	1.3%	12.1%	0.8%	16.1%	-	-
PM-02	0.1%	103.1%	0.4%	105.9%	0.4%	73.3%	14.3%	10.3%	7.0%	3.4%	4.4%	4.4%
PM-03	0.1%	104.2%	0.2%	101.2%	68.0%	12.2%	77.0%	0.6%	66.1%	7.1%	75.6%	1.7%
PM-04	0.1%	103.4%	0.2%	105.5%	0.5%	95.0%	20.2%	42.5%	14.8%	19.9%	2.4%	15.5%
PM-05	0.0%	107.1%	0.0%	106.9%	64.6%	6.6%	70.6%	1.5%	67.5%	0.8%	-	-
PM-07	0.0%	103.3%	1.5%	103.5%	1.9%	95.8%	14.0%	4.7%	4.9%	2.7%	-	-
PM-08	0.0%	108.3%	0.0%	106.5%	2.0%	96.5%	15.5%	15.9%	36.6%	16.6%	-	-
PM-12	0.0%	37.3%	0.2%	103.3%	0.3%	87.3%	48.2%	22.7%	17.0%	10.9%	12.3%	8.9%

BT: Cartridge Break Through

EL: Elution Recovery from Cartridge

