


InertSearch™ for GC

InertCap® Applications

Indoor air by AUTOCAN

Data No. GA011-0000

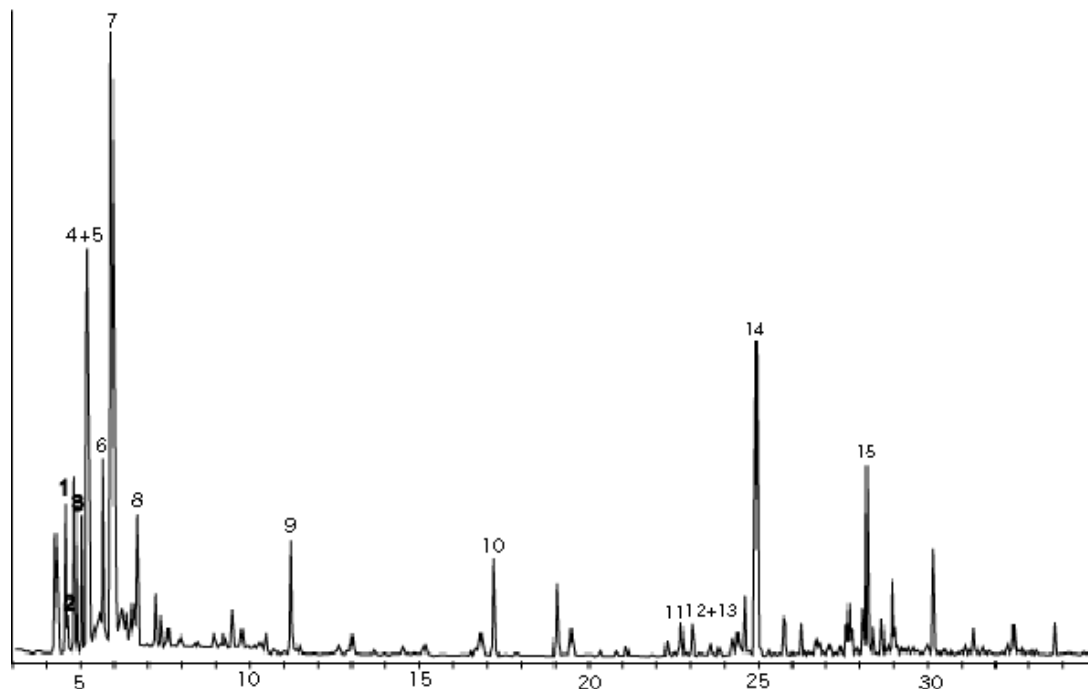

System : GC/MS AUTO Can
Column : InertCap AQUATIC
0.25mmI.D.×60m df=1.0µm
Carrier Gas : He
Injection : AUTO Can/Cryofocus 200mL
Detection : MS
Sample : Indoor air

Solutes : 1. Propane
2. Unknown
3. Butane
4. Methanol
5. Acetaldehyde
6. 2-Methylbutane
7. Ethanol
8. Acetone
9. Ethylacetate
10. Toluene
11. Ethylbenzene
12. m-Xylene
13. p-Xylene
14. α -Pinene
15. Limonene