


## Residual solvents in pharmaceuticals (JP18 Standard stock solution Class 2 B)

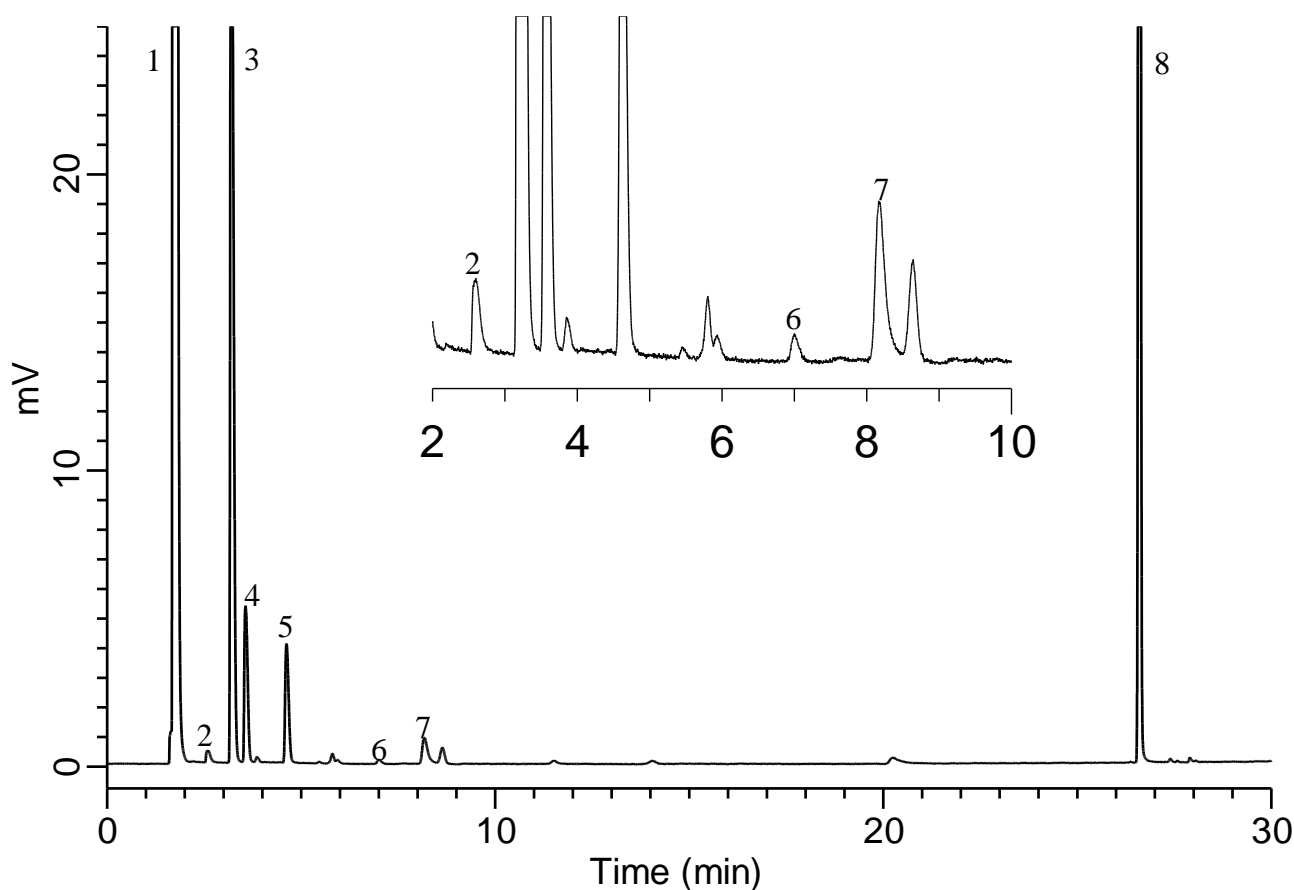

### Conditions

**System** : GC4000 Plus-FID  
**Column** : InertCap WAX (GL Sciences Inc.)  
 0.32 mm I.D. × 30 m df = 0.25 μm  
**Col. Cat. No.** : 1010-67242  
**Col. Temp.** : 50 °C(20 min) – 6 °C/min  
 - 165 °C(20 min)  
**Carrier Gas** : N<sub>2</sub> 2.20 mL/min (35 cm/sec)  
**Injection** : Split 2:1 Manual Injection  
 140 °C  
**Detection** : FID Auto Range  
 250 °C  
**Sample Size** : Head Space Gas 1mL

### Analyte

1. n-Hexane
2. 1,2-Dimethoxyethane
3. Trichloroethylene
4. Chloroform
5. 2-Hexanone(MBK)
6. Nitromethane
7. Pyridine
8. Tetralin