


MonoTrap is a hybrid novel adsorbent that combines a large surface area and the properties of silica gel, activated carbon, and ODS. Utilizing the surface of porous silica, it can be used as a tool that makes use of the large surface area and the adsorbent effect caused by the inclusion of activated carbon for high collection efficiency and for short time and sensitive analysis. In this study, we used MonoTrap DCC18 (with activated carbon) to recover and analyze airborne fatty acids. Activated carbon is effective in collecting polar components such as acid, which can efficiently collect low concentrations of fatty acids. The highly inert WAX-column InertCap Pure-WAX is the optimal column for fragrance components analyses. It is recommended to use this medicine in conjunction with MonoTrap.

Column : InertCap Pure-WAX (Cat.No.1010-68142) 0.25 mm I.D. × 30 m df=0.25 μm

Protocol

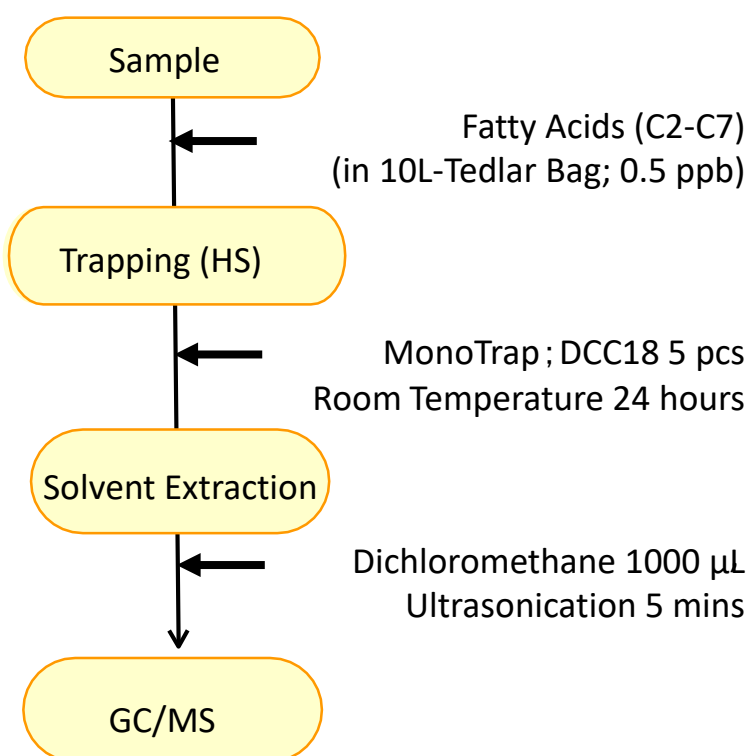

GC Condition

System : SHIMADZU GC-2010, GCMS-QP2010


Column : InertCap Pure-WAX
0.25 mm I.D. × 30 m df=0.25 μm

Column Temp : 50 °C (5 min) → 6 °C/min → 250 °C

Carrier Gas : He 1.15 mL/min


Injection : Split 1:5, 1 μL 250 °C

Detection : MS Scan (*m/z*; 28-450)

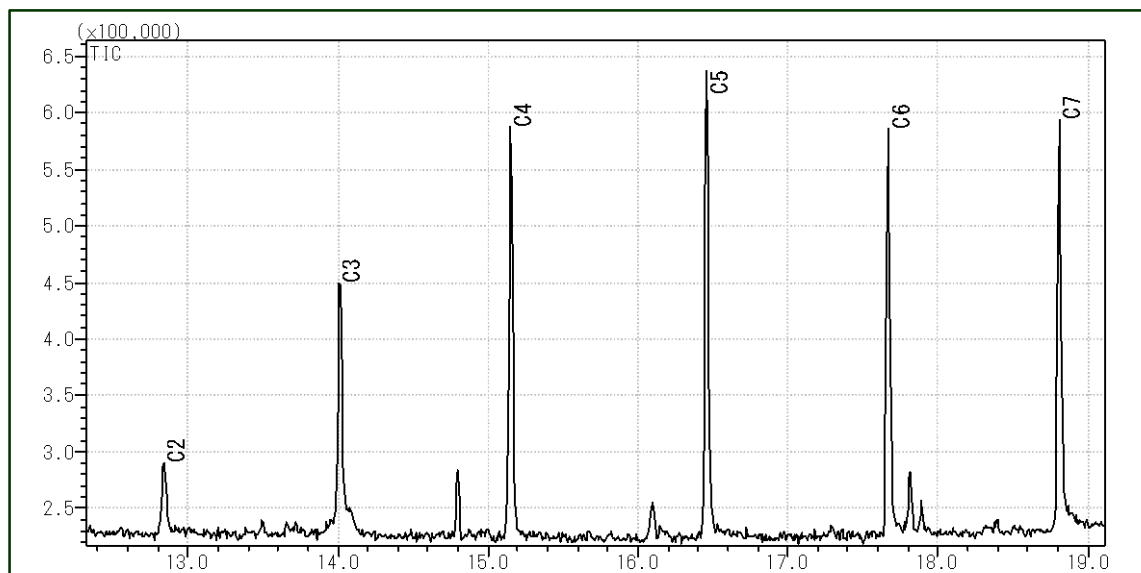

MonoTrap Series

Cat.No	Description	Shape	Size	Active Carbon	Functional Group					
1050-72101	MonoTrap DCC18	Disk	O.D.10 mmx thick 1 mm	Yes	C18					
1050-72201	MonoTrap RCC18	Rod	5mm	Yes	C18					
1050-71101	MonoTrap DSC18	Disk	O.D.10 mmx thick 1 mm	No	1050-71201	MonoTrap RSC18	Rod	5mm	No	C18
1050-71201	MonoTrap RSC18	Rod	5mm	No	C18					

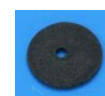

DCC18



RCC18




DSC18



RSC18

MonoTrap Convenient MonoTrapStart-UP-KIT

	Description	Qty.
①	MT Holder	5
②	MT Stand	1
③	MT Extract Cup with Vial (20mL)	5
④	Clean Pin Hole Septum with vial (40mL)	5
⑤	200µL Glass Insert (Flat Bottom)	40
⑥	MonoTrap DCC18	20
⑦	MonoTrap RCC18	20
⑧	MonoTrap DSC18	20
⑨	MonoTrap RSC18	20

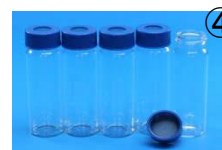

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

GL Sciences, Inc. Japan

22-1 Nishishinjuku 6-Chome
Shinjuku-ku, Tokyo,
163-1130, Japan
Phone: +81-3-5323-6620
Fax: +81-3-5323-6621
Email: world@glsc.co.jp
Web: www.glsciences.com

GL Sciences B.V.

De Sleutel 9
5652 AS Eindhoven
The Netherlands
Phone: +31 (0)40 254 95 31
Email: info@glsciences.eu
Web: www.glsciences.eu

GL Sciences (ShangHai) Ltd.

Tower B, Room 2003,
Far East International Plaza,
NO,317 Xianxia Road,
Changning District.
Shanghai, China P.C. 200051
Phone: +86 (0)21-6278-2272
Email: contact@glsciences.com.cn
Web: www.glsciences.com.cn

GL Sciences, Inc. USA

4733 Torrance Blvd. Suite 255
Torrance, CA 90503
Phone: 310-265-4424
Fax: 310-265-4425
Email: info@glsciencesinc.com
Web: www.glsciencesinc.com

International Distributors

Visit our Website at:
<https://www.glsciences.com/company/distributor.html>

