


Concentration Analysis of VOCs Derived from Expanded Foam Plastic Cushioning Material - Using MonoTrap a Simple Enrichment Tool

MonoTrap DCC18, a simple concentration tool was used to screen and analyze the components that volatilized from expanded foam plastic cushioning material used in a carrier case. The atmosphere in the case was collected using MonoTrap, then extracted with dichloromethane and analyzed by GC/MS. A number of peaks, including alkylbenzene compounds were detected.

Measurement procedure

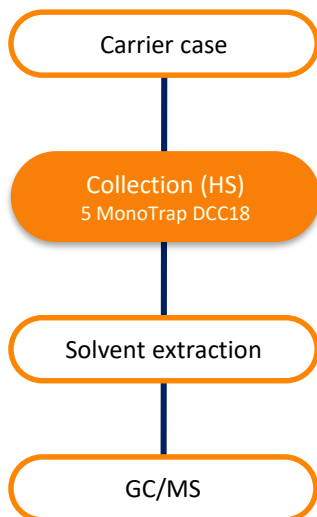

At room temperature for 15 hours




500 μ L of Dichloromethane
5 minutes of ultrasonication

GC/MS Conditions


System	: GC - MS
Column	: InertCap Pure-WAX 0.25 mm I.D. X 30 m, df = 0.25 μ m
Col.Cat. No.	: 1010-68162
Col.Temp.	: 40 °C (5 min) - 4 °C/min - 250 °C
Carrier Gas	: He, 50 kPa
GC Inlet	: 250 °C, Splitless
Sample Size	: 1 μ L
Detection	: MS Scan (m/z 45-450)

GC/MS chromatogram

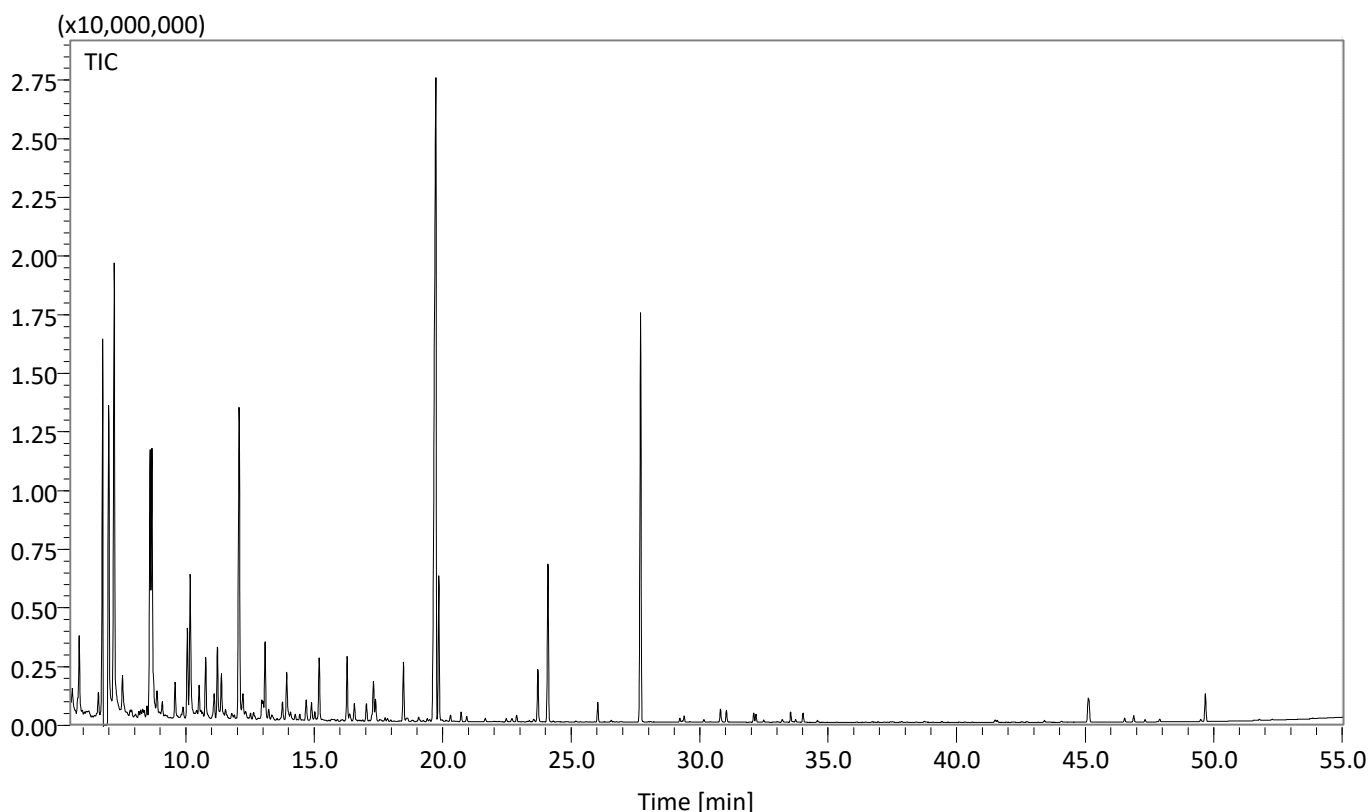

See the following page for the qualitative results of the components obtained by library searching...

Results of library search

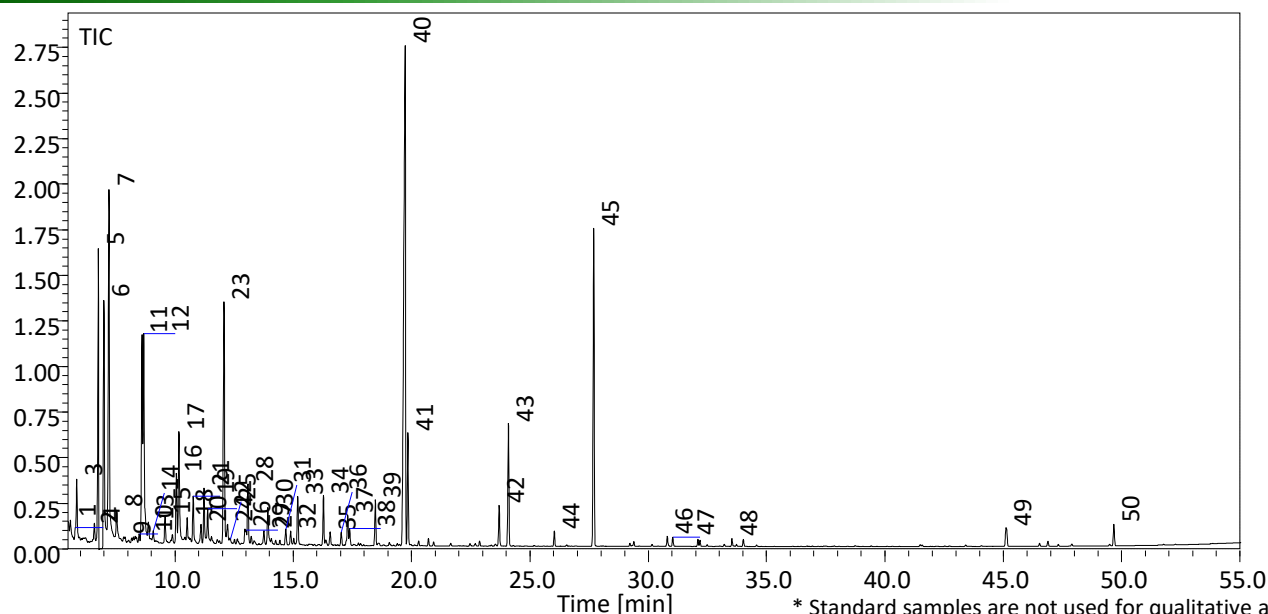

1. Tridecane, 4-cyclohexyl-
2. 1-Propanol, 2-methyl-
3. Propane, 1-(1,1-dimethylethoxy)-2-methyl-
4. Undecane
5. Ethylbenzene
6. Xylene
7. Xylene
8. 1-Butanol
9. 1-Undecene
10. 2,3-Dimethyldecane
11. Xylene
12. 4-Heptanone, 2,6-dimethyl-
13. Undecane, 4-cyclohexyl-
14. Trans-Decalin, 2-methyl-
15. Benzene, propyl-
16. Benzene, 1-ethyl-2-methyl-
17. Benzene, 1-ethyl-3-methyl-
18. 1-Methoxy-2-propyl acetate
19. Benzene, 1,3,5-trimethyl-
20. 2-Heptanone, 4,6-dimethyl-
21. Styrene
22. Benzene, 1-ethyl-2-methyl-
23. Benzene, 1,3,5-trimethyl-
24. Propanoic acid, 2-hydroxy-, 2-methylpropyl ester
25. 9-Eicosyne

26. Benzene, 1-methyl-2-(1-methylethyl)-
27. Benzene, 1-methyl-2-propyl-
28. 2-Ethoxyethyl acetate
29. Benzene, 1-ethyl-3,5-dimethyl-
30. Benzene, 1,2,3-trimethyl-
31. Benzene, 1-methyl-2-(1-methylethyl)-
32. Benzene, 1-methyl-3-(1-methylethyl)-
33. Benzene, 2-ethyl-1,4-dimethyl-
34. Acetic acid, 2-ethylhexyl ester
35. Ethanol, 2-butoxy-
36. Benzene, 1,2,4,5-tetramethyl-
37. Benzene, (1-methoxy-1-methylethyl)-
38. Benzene, 1,2,3,5-tetramethyl-
39. 2-Ethylhexyl chloroformate
40. 2-Propenoic acid, 2-ethylhexyl ester
41. 1-Hexanol, 2-ethyl-
42. Tetramethylbutanedinitrile
43. Acetophenone
44. Pentanedioic acid, dimethyl ester
45. Benzenemethanol, -dimethyl-
46. Propanoic acid, 2-methyl-, 3-hydroxy-2,4,4-trimethylpentyl ester
47. 2,2,4-Trimethyl-1,3-pentanediol diisobutyrate
48. 2H-Pyran-3(4H)-one, 6-ethenyldihydro-2,2,6-trimethyl-
49. Tetramethyl-1,4-benzenedimethanol
50. Dibutyl phthalate

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

GL Sciences, Inc. Japan

22-1 Nishishinjuku 6-Chome
Shinjuku-ku, Tokyo,
163-1130, Japan
Phone: +81-3-5323-6620
Fax: +81-3-5323-6621
Email: world@glsc.co.jp
Web: www.glsciences.com

GL Sciences B.V.

De Sleutel 9
5652 AS Eindhoven
The Netherlands
Phone: +31 (0)40 254 95 31
Email: info@glsciences.eu
Web: www.glsciences.eu

GL Sciences, Inc. USA

4733 Torrance Blvd. Suite 255
Torrance, CA 90503
Phone: 310-265-4424
Fax: 310-265-4425
Email: info@glsciencesinc.com
Web: www.glsciencesinc.com

GL Sciences (ShangHai) Ltd.

Tower B, Room 2003,
Far East International Plaza,
NO,317 Xianxia Road,
Changning District.
Shanghai, China P.C. 200032
Phone: +86 (0)21-6278-2272
Email: contact@glsciences.com.cn
Web: www.glsciences.com.cn

International Distributors

Visit our Website at www.glsciences.com/distributors