


## Analysis of Engine oil

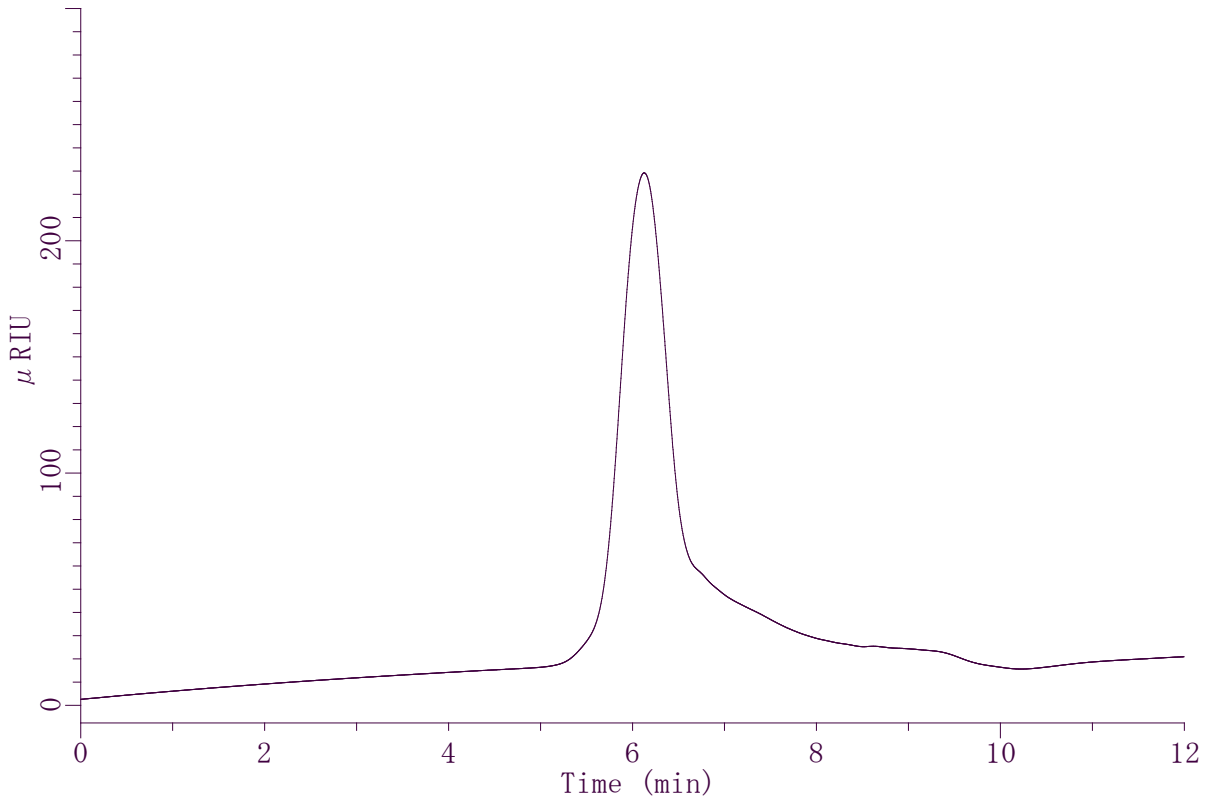

### Conditions

**System** : GL7400 HPLC system  
**Column** : Inertsil Diol (5 $\mu$ m, 250 x 7.6 mmI.D.)  
**Column Cat. No.** : 5020-05666  
**Eluent** : THF  
**Flow Rate** : 1.0 mL/min  
**Col. Temp.** : 40°C  
**Detection** : RI (GL-7454 RI Detector) (Cell Temp. 35°C)  
**Injection Vol.** : 10 $\mu$ L  
**Sample** : Engine oil