
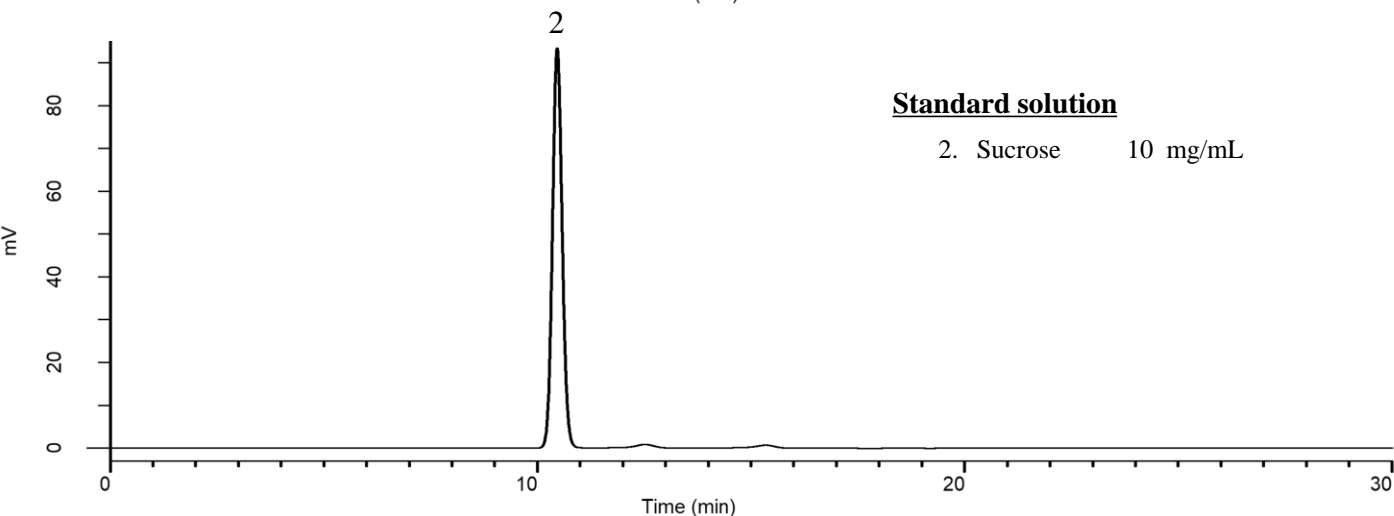


Analysis of Sucrose

(Under the condition of the PDG (Pharmacopeial Discussion Group) Stage 2 Drafts)

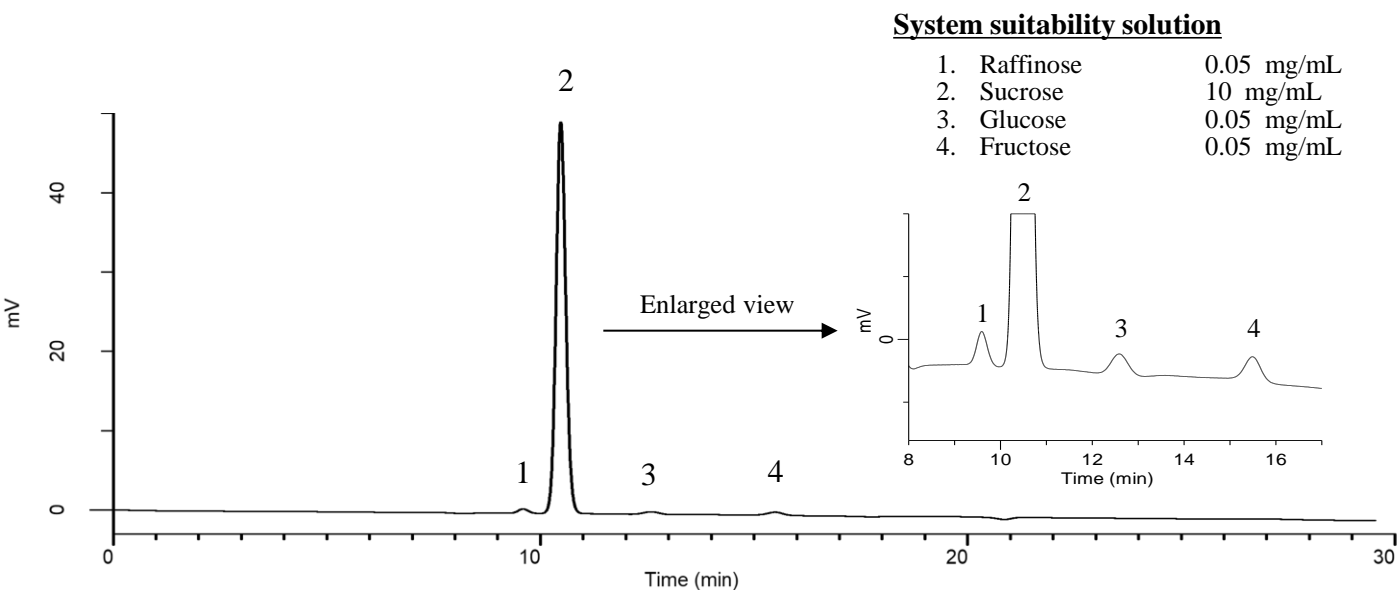

Conditions

System : Chromaster
Column : InertSphere Sugar-2
 (9 μ m, 300 x 7.8 mm I.D.)
Column Cat. No. : 5020-11000
Eluent : H₂O
Flow Rate : 0.5 mL/min
Col. Temp. : 80 °C
Detection : RI (5450 RI Detector)
Injection Vol. : 10 μ L
Sample : Standard

Analyte:

1. Raffinose : Relative retention time (1 to 2) : 0.92 (about 0.9)
 2. Sucrose
 3. Glucose : Relative retention time (3 to 2) : 1.19 (about 1.2)
 4. Fructose : Relative retention time (4 to 2) : 1.48 (about 1.6)

System suitability solution

Peak-to-valley ratio (1, 2) : 31.07 (\geq 2.5)
 Resolution (2, 3) : 3.86 (\geq 1.5)

Standard solution

RSD of the Peak area of 2 (%) (n=5) : 0.11 (\leq 0.73)