

Development of polar lipid profiling method by supercritical fluid chromatography/mass spectrometry

Data No. LL007-0000

The poster was presented at The 59th Annual Conference on Mass Spectrometry, Osaka, Japan, 2011.

Provided by Dr. Takeshi Bamba, an associate professor, Department of Biotechnology, Graduate School of Engineering Osaka University, 2-1 Yamadaoka, Suita, Osaka 565-0871, Japan

Author name : Jae Won Lee; Takashi Yamamoto; Takato Uchikata; Atsuki Matsubara; Eiichiro Fukusaki; Takeshi Bamba
Data source : Poster
Year : 2011

Conditions

Column : Inertsil ODS-EP (5 μm, 250 x 4.6 mm I.D.)
Column Cat. No. : 5020-02646
Mobile Phase : Carbon dioxide (CO₂, 99.9% grade)
Modifier : Methanol with 0.1% (w/w) HCOONH₄
Flow rate : 3 mL/min
Oven temperature : 37 °C
Back pressure : 10 MPa

Sample : Polar lipids
Analyte : table below

Detection : SFC/MS/MS
Ionization method : Electrospray ionization
Ion mode : positive
Capillary voltage : 3.00 kV
Desolvation temperature : 450 °C
Desolvation gas flow : 800 L/hr
Cone gas flow : 60 L/hr
Collision gas flow : 13.2 mL/hr
MS collision energy : 6 V
Source temperature : 150 °C
Extractor voltage : 3 V

Optimized MS/MS method by TMS derivatization

Polar lipids	SRM transitions	CV	CE	Number of adducted TMS
Phosphatidylglycerol (PG)	[M+H] ⁺ > [M-315]	20	30	2
Phosphatidic acid (PA)	[M+H] ⁺ > [M-169]	25	25	1
Phosphatidylinositol (PI)	[M+H] ⁺ > [M-619]	35	30	5
Lysophosphatidylcholine (LPC)	[M+H] ⁺ > 184	45	35	1
Lysophosphatidylethanolamine (LPE)	[M+H] ⁺ > [M-140]	30	15	1
Lysophosphatidylglycerol (LPG)	[M+H] ⁺ > [M-315]	30	20	3
Lysophosphatidic acid (LPA)	[M+H] ⁺ > [M-169]	30	15	2
Lysophosphatidylinositol (LPI)	[M+H] ⁺ > [M-547]	45	25	5
Sphingomyeline (SM)	[M+H] ⁺ > 184	45	30	1
Sphingosine-1-phosphate (SIP)	[M+H] ⁺ > 264	25	20	2

CV: Cone voltage, CE: MS/MS collision energy

Profiling of 10 polar lipid standards

