
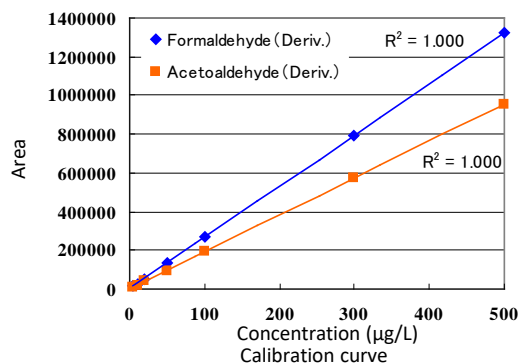


Aldehydes are widely recognized as hazardous pollutants in both the atmosphere and building interiors. This report introduces the methods of analysis for formaldehyde and acetaldehyde, which are two of the most commonly analyzed aldehydes. A High Performance Liquid Chromatography (HPLC) was used as an analyzer.

Aldehydes were collected using the "sampling bag method," which is used mainly for VOC measurement in automotive interior materials.

As a pretreatment, the sample is placed in a Tedlar bag and heated for a period of time to release the aldehydes. The gas is then suctioned out with a sampling pump and collected on a DNPH cartridge. Acetonitrile is passed through the DNPH collection cartridges to elute the aldehydes. The recovered eluent is analyzed by HPLC to detect the aldehydes.

## Example: Measurement of standard solution

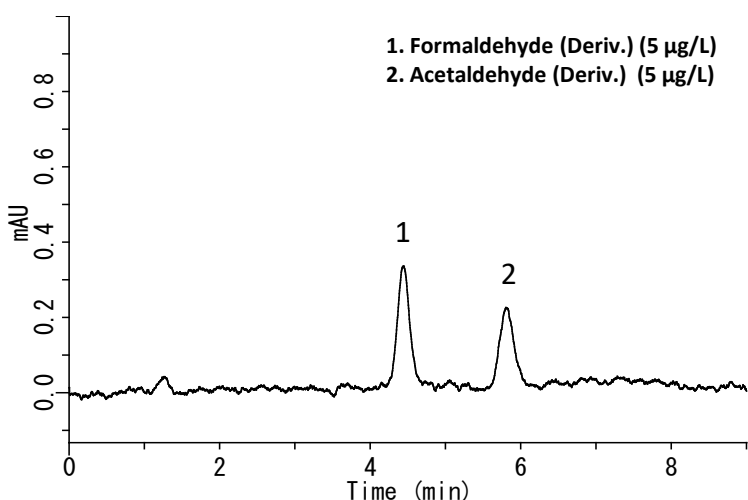

### HPLC conditions

<b>Column</b>	: Inertsil ODS-80 A (5 µm, 150 x 4.6 mm I.D.)
<b>Eluent</b>	: A) CH <sub>3</sub> CN B) H <sub>2</sub> O A/B = 55/45, v/v (gradient mixer)
<b>Flow rate</b>	: 1.0 mL/min
<b>Column temperature</b>	: 40 °C
<b>Detected</b>	: UV 360 nm
<b>Injection volume</b>	: 20 µL

## Measurement example

### Sample pretreatment example

#### Sample

- Place the sample piece in a Tedlar-bag.
- Include a fixed amount of nitrogen gas

#### Heating

- At -60 °C to 65 °C and heated for 2 h.


#### Collecting

- Collect in DNPH cartridges

#### Collection


- Acetonitrile 5 mL

#### Measurement

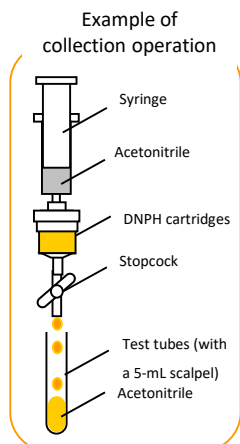


Measurements were made using the sampling bag-method and HPLC.

### Resin

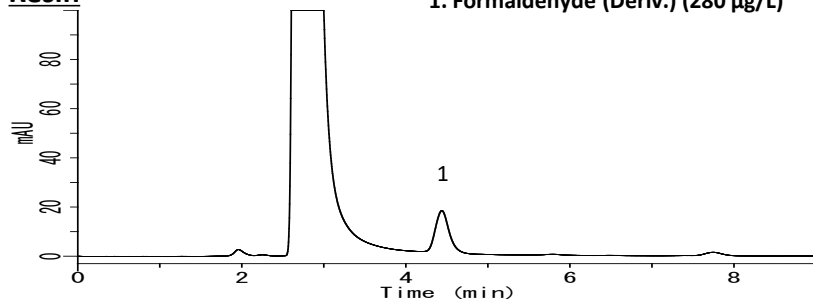

### Car interior materials

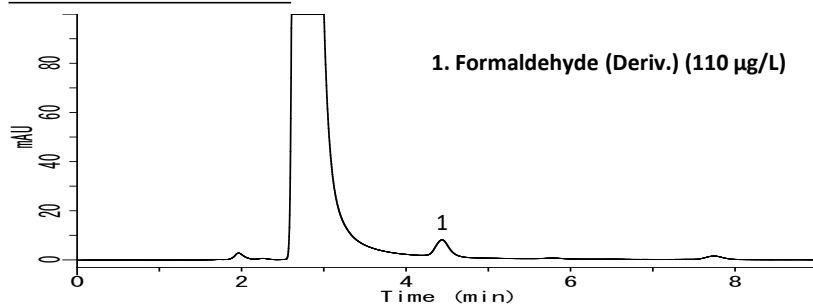

## Overview of the sampling system

This section details the system for recovering aldehyde gases released from materials using the "Sampling Bag Method."

### Aldehyde sampler: sampler bag method-compatible sampler set

#### <Features of the sampling bag method collection system>

- Two samples can be heated and collected simultaneously.
- It can also be used for VOC collection.
- A nitrogen gas supply unit is included

#### <DNPH cartridges: GL-Pak mini AERO DNPH>


- Carbonyl compounds are collected using silica gel impregnated with 2,4-dinitrophenylhydrazine (DNPH).
- Due to the use of spherical silica gel, this cartridge has excellent air permeability, high collection efficiency, and low blank.

#### <Sampling pumps: SP208-Dual 1000>

- Two lines can be used simultaneously.
- A data log can be recorded and is useful for quality control.
- Even if the suction resistance fluctuates, the set flow rate can be maintained and suction is made at a stable rate.

#### <Tedlar bag>

- Plastic bags are available for a wide range of applications, from inorganic gases to organic solvents.

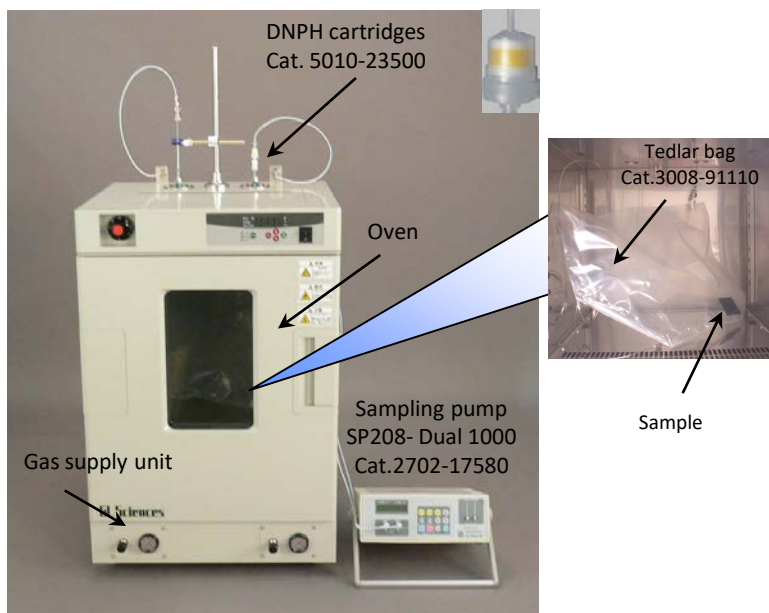

Sampling bag method collection system

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

#### **GL Sciences, Inc. Japan**

22-1 Nishishinjuku 6-Chome  
Shinjuku-ku, Tokyo,  
163-1130, Japan  
Phone: +81-3-5323-6620  
Fax: +81-3-5323-6621  
Email: [world@glsciences.co.jp](mailto:world@glsciences.co.jp)  
Web: [www.glsciences.com](http://www.glsciences.com)

#### **GL Sciences B.V.**

De Sleutel 9  
5652 AS Eindhoven  
The Netherlands  
Phone: +31 (0)40 254 95 31  
Email: [info@glsciences.eu](mailto:info@glsciences.eu)  
Web: [www.glsciences.eu](http://www.glsciences.eu)

#### **GL Sciences (Shanghai) Ltd.**

Tower B, Room 2003,  
Far East International Plaza,  
NO,317 Xianxia Road,  
Changning District.  
Shanghai, China P.C. 200032  
Phone: +86 (0)21-6278-2272  
Email: [contact@glsciences.com.cn](mailto:contact@glsciences.com.cn)  
Web: [www.glsciences.com.cn](http://www.glsciences.com.cn)

#### **GL Sciences, Inc. USA**

4733 Torrance Blvd. Suite 255  
Torrance, CA 90503  
Phone: 310-265-4424  
Fax: 310-265-4425  
Email: [info@glsciencesinc.com](mailto:info@glsciencesinc.com)  
Web: [www.glsciencesinc.com](http://www.glsciencesinc.com)

#### **International Distributors**

Visit our Website at [www.glsciences.com/distributors](http://www.glsciences.com/distributors)