


$\alpha$ -Carotene and  $\beta$ -carotene are a type of carotenoid and widely known as precursors to vitamin A in human body. They also have antioxidant activity.

The two carotenes are very similar in chemical structure. Since the difference is only in the position of a carbon-carbon double bond, it is difficult to separate them with widely-used monomeric ODS columns. To solve this problem, Inertsil ODS-P, which is a polymeric ODS column and superior in steric recognition, was used as a separation column in this note.



(K. Suzuki)

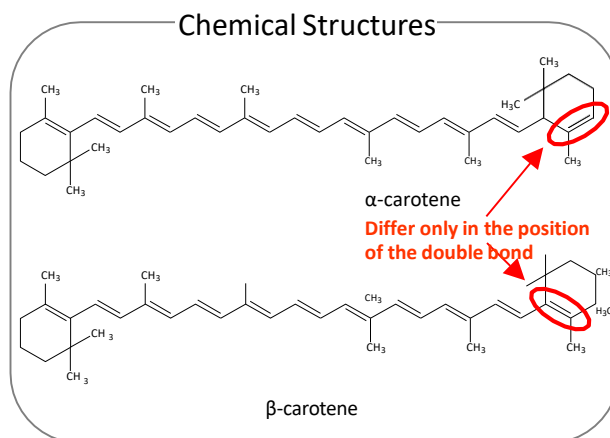

### Chromatograms obtained from standard solution

When a commercially available monomeric column was used,  $\alpha$ -carotene and  $\beta$ -carotene were not separated within 20 minutes. However, Inertsil ODS-P provided good separation of the two carotenes.


1.  $\alpha$ -Carotene (2.0 mg/L)
2.  $\beta$ -Carotene (2.0 mg/L)

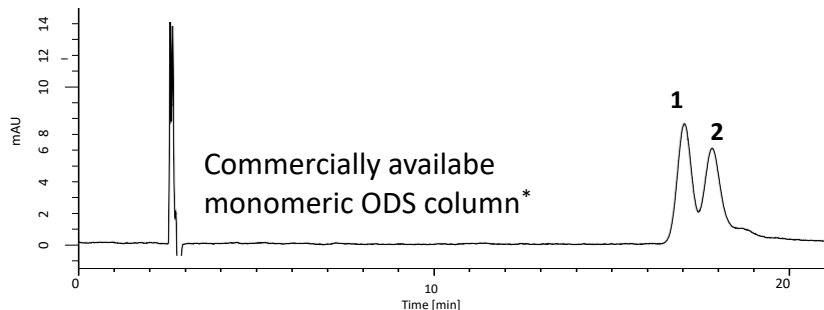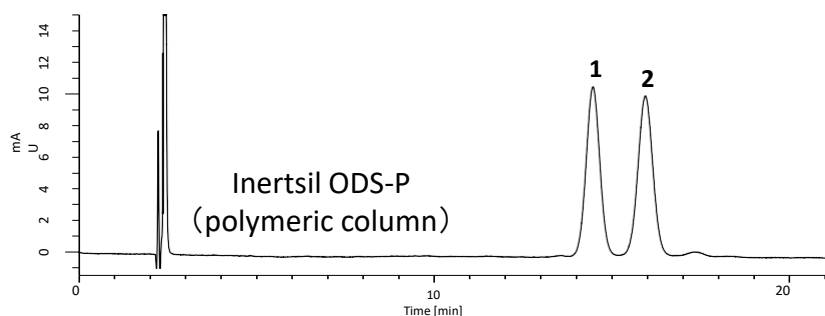

\* The mobile phase composition was adjusted to A/B = 45/55.

### Conditions

- Column** : Inertsil ODS-P  
(5  $\mu$ m, 250 x 4.6 mm I.D.)  
Cat.No. 5020-02002
- Eluent** : A) CH<sub>3</sub>OH  
B) C<sub>2</sub>H<sub>5</sub>OH  
A/B = 90/10, v/v  
(gradient mixer)
- Flow rate** : 1.0 mL/min
- Col. Temp.** : 40 °C
- Detection** : PDA 455 nm
- Inj. Vol.** : 20  $\mu$ L

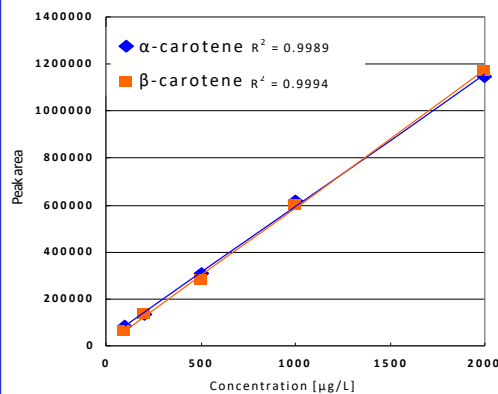

Calibration curves obtained with  
Inertsil ODS-P column

## Analyses of food samples

### Sample pretreatment 1

#### Vegetables

- 0.5 g
- 3 % Pyrogallol ethanol solution 10 mL
- 60 % KOH aqueous solution 1 mL

#### Saponification

- Heat at 70 °C for 30 min
- Place in an ice-water bath to cool down
- 1 % NaCl aqueous solution 22.5 mL
- Hexane-ethyl acetate = 9/1 (v/v) 15 mL
- Shaking for 5 min

#### Liquid-liquid extraction

- Organic layer
- Aqueous layer
- Repeat procedure \*A twice

#### Concentration

- Evaporate under reduced pressure
- Dissolve in 2 mL of ethanol

#### HPLC

### Sample pretreatment 2

#### Carrot juice

- 2 g
- 3 % Pyrogallol ethanol solution 20 mL
- Sodium sulfate anhydrous 10 g
- Shaking for 5 min

#### Centrifuge

- Supernatant
- Residue
- Repeat procedure \*B twice
- 1 % NaCl aqueous solution 23 mL
- 2-Propanol 6 mL
- Hexane-ethyl acetate = 9/1 (v/v) 15 mL
- Shaking for 5 min



#### Liquid-liquid extraction

- Organic layer
- Aqueous layer
- Repeat procedure \*A twice

#### Concentration

- Evaporate under reduced pressure
- Dissolve in 10 mL of ethanol

#### HPLC

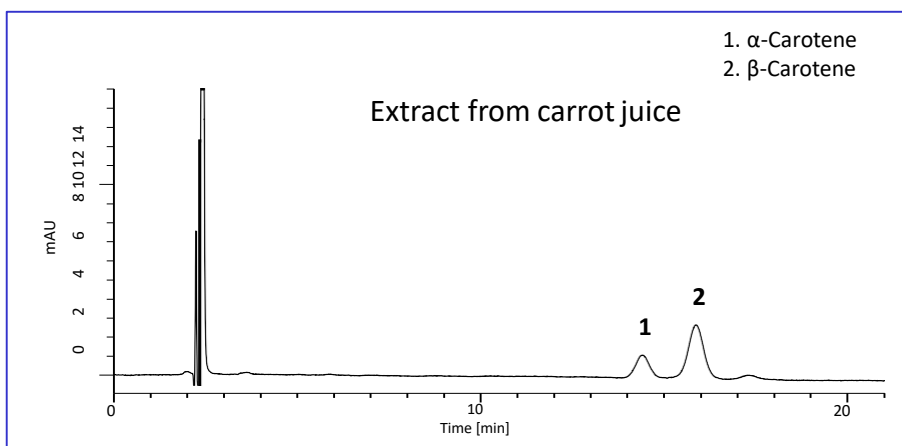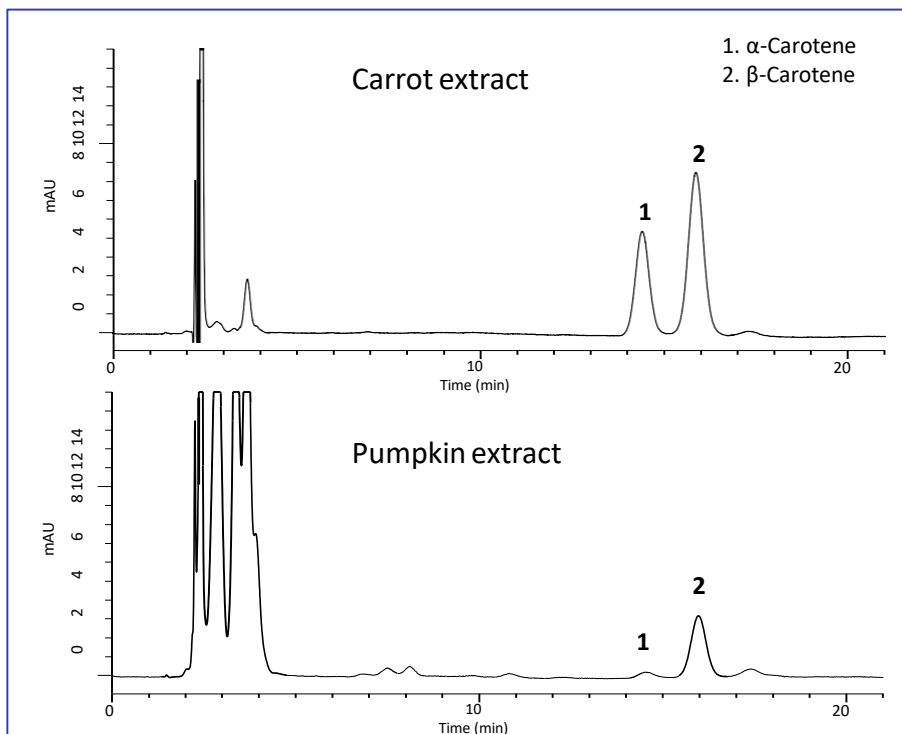

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

#### GL Sciences, Inc. Japan

22-1 Nishishinjuku 6-Chome  
Shinjuku-ku, Tokyo,  
163-1130, Japan  
Phone: +81-3-5323-6620  
Fax: +81-3-5323-6621  
Email: [world@glsc.co.jp](mailto:world@glsc.co.jp)  
Web: [www.glsciences.com](http://www.glsciences.com)

#### GL Sciences B.V.

De Sleutel 9  
5652 AS Eindhoven  
The Netherlands  
Phone: +31 (0)40 254 95 31  
Email: [info@glsciences.eu](mailto:info@glsciences.eu)  
Web: [www.glsciences.eu](http://www.glsciences.eu)

#### GL Sciences (ShangHai) Ltd.

Tower B, Room 2003,  
Far East International Plaza,  
NO,317 Xianxia Road,  
Changning District.  
Shanghai, China P.C. 200032  
Phone: +86 (0)21-6278-2272  
Email: [contact@glsciences.com.cn](mailto:contact@glsciences.com.cn)  
Web: [www.glsciences.com.cn](http://www.glsciences.com.cn)

#### GL Sciences, Inc. USA

4733 Torrance Blvd. Suite 255  
Torrance, CA 90503  
Phone: 310-265-4424  
Fax: 310-265-4425  
Email: [info@glsciencesinc.com](mailto:info@glsciencesinc.com)  
Web: [www.glsciencesinc.com](http://www.glsciencesinc.com)

#### International Distributors

Visit our Website at:  
<https://www.glsciences.com/company/distributor.html>